Mikroskop · Microscope ·

Science TFM-201/301 + Science Infinity

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EN Instruction manual

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Contents

1 Imprint	16
2 Validity information	16
3 About this Instruction Manual.....	16
4 Intended Use	16
5 General safety instructions	17
6 Product images.....	19
7 Scope of delivery	19
8 Parts overview	19
9 Site location	20
10 Installing the USB power adapter	20
11 Setting up power supply.....	21
12 Turning on/off the device.....	21
13 Adjusting the lighting.....	21
14 Inserting slides	22
15 Observation.....	22
16 Adjusting the eye distance	23
17 Diopter adjustment.....	23
18 Using the iris diaphragm	23
19 Using the Köhler illumination.....	23
20 Oil immersion.....	24
21 Using the trinocular extension	24
22 Cleaning and maintenance	24
23 Disposal.....	25
24 Technical data.....	25
25 Warranty	26
26 EC Declaration of Conformity	26

1 Imprint

Bresser GmbH

Gutenbergstr. 2

46414 Rhede

Germany

<http://www.bresser.de>

For information regarding liability claims or service requests, please refer to chapters "Warranty" and "Service" in this documentation. We ask for your understanding, that any requests or submissions that have been addressed to the manufacturer's address can not be processed.

Errors reserved - technical specifications subject to change.

Copyright: 2017Bresser GmbH

All rights reserved.

It is not allowed to reproduce this documentation or parts of it in any kind (e.g. photocopying, printing, etc) without written a permission of the manufacturer. It is also prohibited to use or transmit this documentation with electronic systems (e.g. image file, website, etc).

The descriptions and brand names of the corresponding firms used in this documentation are generally protected under trademark, brand and/or patent law for Germany, for the European Union and/or other countries.

2 Validity information

This documentation is valid for the products with the article numbers listed below:

5750900

Manual version:v032017a

Manual description:Manual_Science-TFM-Infinity-Series_de-en_BRESSER_v032017a

With any service inquiries, please state these information.

3 About this Instruction Manual

NOTICE

These operating instructions are to be considered a component of the device.

Please read the safety instructions and the operating instructions carefully before use.

Keep these instructions for renewed use at a later date. When the device is sold or given to someone else, the instruction manual must be provided to the new owner/user of the product.

4 Intended Use

- This device is not intended for use by individuals (including children) with limited physical, sensory or mental capabilities or those lacking in experience and/or knowledge, unless they are supervised by an individual responsible for their safety or have received instructions from them regarding the use of the device.
- This product is intended only for private use.
- It was developed for the magnified display of things in nature.
- The device is intended only for indoor use.

5 General safety instructions

DANGER

Danger of an electric shock!

This device contains electronic components that operate via a power source (power supply and/or batteries). In case of any improper use of this device, there is a risk of an electric shock. An electric shock can cause severe injury or even death. Therefore please read the safety instructions below to avoid an explosion.

- Disconnect the device from the power supply by pulling the power plug when it is not used or in case of longer interruption of operation and before starting any work on maintenance and cleaning.
- Position your device so that it can be disconnected from the power supply at any time. The wall socket you use should be located near the device and easily accessible, since the plug on the power cable serves as a disconnecting device for the power supply.
- Always pull on the plug to separate the device from the power supply. Never pull on the cable.
- Before operating, check the device, cables and connections for damage.
- Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.
- Only use the device in complete dry environment and do not touch it with wet or moist parts of your body.

DANGER

Danger of choking!

In case of any improper use of this device, there is a risk of choking, especially for children. Therefore please read the safety instructions below.

- Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard!

DANGER

Danger of explosion!

In case of any improper use of this device, there is a risk of an explosion. Therefore please read the safety instructions below to avoid an explosion.

- Do not expose the device to high temperatures. Use only the supplied power adapter. Do not short-circuit the device or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.

CAUTION

Danger of injury!

This device contains components and/or accessories that can cause minor to severe injuries in case of any improper use. Therefore please read the safety instructions below to avoid any bodily injury.

- Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.

-
- Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after using the chemicals. In the event that the chemicals come into contact with your eyes or mouth, rinse thoroughly with water. If you are in pain after exposure, contact a doctor immediately and take the substances with which you came into contact with you.

CAUTION

Fire hazard!

In case of any improper use of this device, there is a risk of fire. Therefore please read the safety instructions below to avoid the initiation of burning.

-
- Never cover the ventilation slots or cooling fins of the device while using it or as long as it has not sufficiently cooled down!

NOTICE

Risk of property damage!

In case of any improper use of this device and/or its accessories, there is a risk of property damage. Therefore only use the device according to the safety instructions below.

-
- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
 - Do not expose this device to higher temperatures and protect it from water and high humidity.
 - Protect the device from severe shocks!
 - For this device only use accessories and spare parts that comply with the technical information.

6 Product images

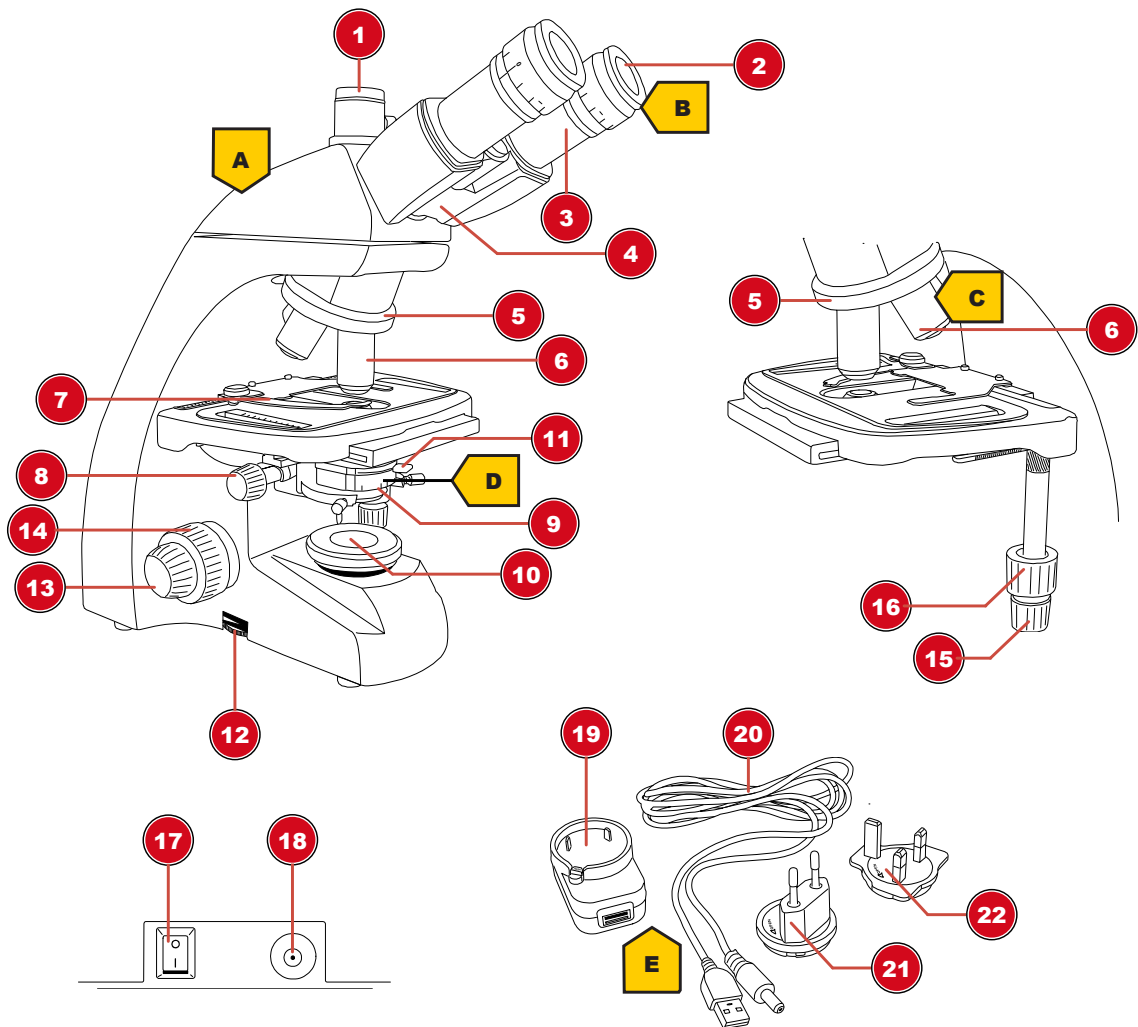
Illustration 2: Product images

7 Scope of delivery

Microscope (A); 10x WF Eyepiece (B); 4 Objectives (pre-assembled) (C); Condenser (pre-assembled) (D); Power adapter with plug and cable (E)

8 Parts overview

1	Trinocular head* / photo adapter*
2	Eyepiece
3	Eyepiece holder
4	Microscope head
5	Objective nosepiece
6	Objective
7	Specimen clamp
8	Condenser height adjustment
9	Condenser

10	Illumination
11	Adjusting lever for iris diaphragm
12	Rotary knob for brightness adjustment
13	Fine focusing
14	Coarse focusing
15	Cross table lateral adjustment (X-axis)
16	Cross table longitudinal adjustment (Y-axis)
17	On/Off switch
18	DC power port
19	Universal USB power adapter
20	USB power cable with coaxial/barrel connector
21	Power adapter attachment (EU)
22	Power adapter attachment (UK)

*not for model TFM-201

9 Site location

Choose a suitable location before assembly and commissioning.

Place the device on a stable and flat surface, free from vibrations.

10 Installing the USB power adapter

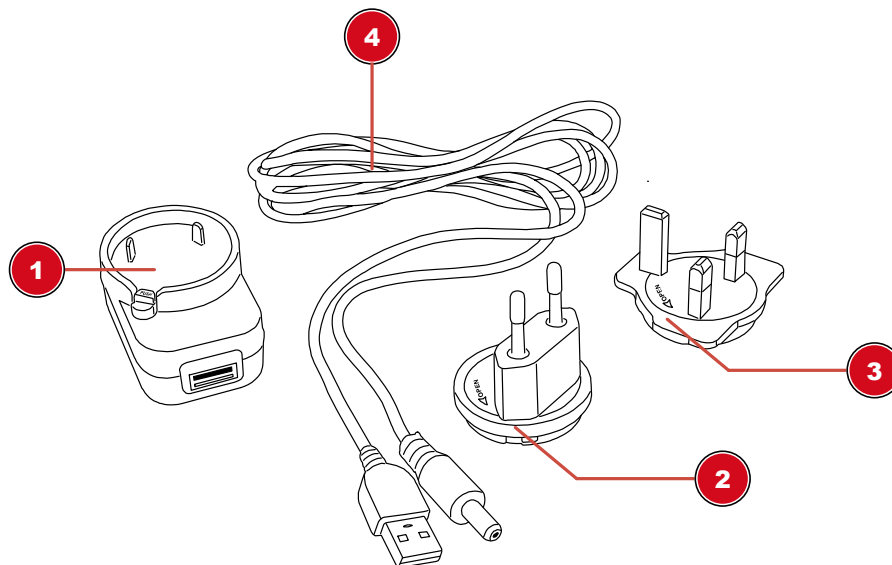
Illustration 3: USB power adapter with exchangeable attachment

1 USB power adapter	2 Adapter attachment (EU)
3 Adapter attachment (UK)	4 USB power cable with coaxial/barrel connector

1. Place suitable adapter attachment with its notches on the pins of the USB power adapter.
2. Turn adapter attachment in LOCK direction until it is locked.
3. Put the USB plug of the power cable into the USB port of the power adapter.
4. To remove the adapter attachment, push the PUSH button and turn the attachment in OPEN direction.

NOTICE

Using the USB power cable with a computer

Alternatively the device can be powered by a computer via the included USB power cable: Put the USB plug into a free USB port of the computer.

11 Setting up power supply

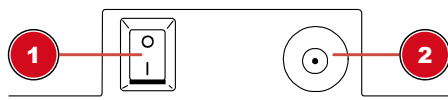
Illustration 4: DC connection socket for coaxial/barrel connector

1 On/Off switch

2 DC connection socket

1. Put the DC connector of the power cable into the DC connection socket of the device.
2. Put the mains plug of the power cable into the power outlet.

12 Turning on/off the device

1. Move On/Off switch to position 'I' to turn on the device.
2. Move On/Off switch to 'O' position to turn off the device.

13 Adjusting the lighting

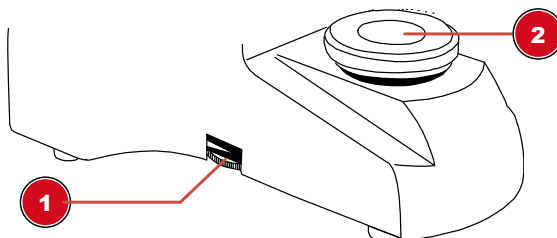
Illustration 5: Lighting unit

1 Rotary knob for brightness adjustment

2 Illumination

This device provides an illumination with a permanently installed LED light source.

NOTICE

Do not operation the device on highest brightness over a longer time!

This will reduce the lifetime of the LED.

To control the brightness proceed as follows:

1. Turn on the device.
2. Turn/Move the dimmer control to adjust the brightness of the light source.

NOTICE

Illuminating the object optimally

Adjust the brightness of the light source so that the illumination works at its optimum. Depending on light influences from outside and the object being observed, the illumination needs regular adjustments.

14 Inserting slides

1. Move the object plate downwards completely by turning the focus wheel.
2. Press the object clamping outwards.
3. Insert the slide straight into the fixture's angle.
4. Let the object clamping fall back carefully until the slide is fixed in the fixture.

15 Observation

NOTICE

Using eyepieces and objectives

When using a specific eyepiece or objective, you will get the individual magnification of an object being observed. The magnification is imprinted on the eyepiece and objective with a large number (e.g. 10x for 10-fold magnification, 4 for 4-fold magnification).

1. Move the object plate downwards completely by turning the focus wheel.
2. Insert a slide and place it in the middle under the objective.
3. Set an objective with lowest magnification below the object plate.
4. Insert an eyepiece with lowest magnification in the eyepiece holder.
5. Turn on the device and the illumination.
6. Look through the eyepiece and move the object plate upwards by turning the focus wheel until a sharp image is visible.
7. If necessary, adjust the image sharpness by turning the fine focus adjustment.

NOTICE

Avoid damages to the specimen!

Only turn the focus wheel until it stops to avoid damages to the specimen throughout collision with the objective.

16 Adjusting the eye distance

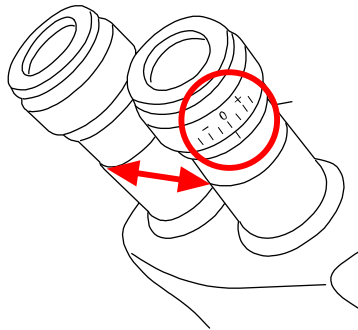
Illustration 6: Distance between the eyepieces and diopter adjustment

1. With the eyes look through both eyepieces.
2. Change the distance between the eyepieces by pushing them together or pulling apart until a circular image of the observed object is visible.

17 Diopter adjustment

1. Turn the diopter adjustment ring to position '0'.
2. With the left eye, look through the left eyepiece and turn the focusing wheel until a sharp image of the observed object is visible.
3. With the right eye, look through the right eyepiece and turn the diopter adjustment ring until a sharp image of the observed object is visible too.

The diopter adjustment can also be done in the opposite way on the left eyepiece.

18 Using the iris diaphragm

NOTICE

The iris diaphragm is used to increase the depth of focus.

With an open aperture more light is getting through but the depth of focus is less. With partially closed aperture the depth of focus is higher but it is necessary to increase the brightness of the illumination then.

It is recommended not to close the iris diaphragm too much when working with higher magnifications. Otherwise the image resolution will decrease.

Move the lever for the iris diaphragm to right or left to change the aperture setting.

19 Using the Köhler illumination

Proceed as follows to adjust the Köhler illumination for an optimal illumination of the field of view:

1. Focus on the object with the 10x objective.
2. Close the condensor and the light field aperture. For microscopes with additional lens, fold it out.
3. Move the condensor up and down until the edge of the light field aperture is in focus.
4. If necessary, use the adjustment screws to move the condensor until the light field aperture is centered in the field of view.
5. Open the light field aperture as far as it is not visible in the field of view anymore.

20 Oil immersion

The oil immersion can be done with a suitable 100x objective only. For the purpose proceed as follows:

1. Focus on the object with the 40x objective.
2. Turn the 40x objective sideways.
3. Drop a drop of immersion oil onto the cover glass of the preparation.
4. Retract the 100x objective so that it is immersed in the oil.
5. Refocus the image.
6. Clean the 100x objective after each use.

21 Using the trinocular extension

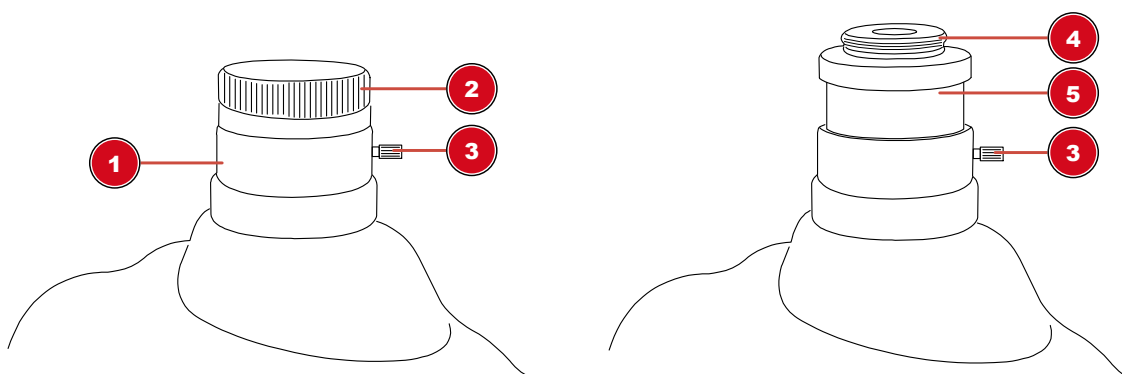
Illustration 7: Trinocular extension with pre-assembled photo adapter

1 Trinocular extension	2 Protection cap
3 Fixing screw	4 Photo adapter thread
5 Photo adapter	

The The trinocular extension of the microscope was specially designed for C-mount threads. For the camera mounting proceed as follows:

1. Remove the dust protection cap from the trinocular head.
2. Attach the camera* with its C-mount thread to the photo adapter thread and turn it on.
3. Loosen the clamping screw of the trinocular extension to move the photo adapter with the camera* within the extension tube. In this way the image sharpness and the position of the camera can be adjusted. Most camera models can be adjusted parfocal to the eyepieces when doing so.
4. Re-tighten the screw to fix the camera in the desired position.

*not included

22 Cleaning and maintenance

- Protect the device against dust and moisture.
- Avoid putting fingerprints and equal contaminations on any optical surfaces.
- Make sure the residual moisture is fully eliminated before use.
- Disconnect the device from the power supply by pulling the power plug when it is not used or in case of longer interruption of operation and before starting any work on maintenance and cleaning.
- Only use a dry cloth to clean the exterior of the device. To avoid damaging the electronics, do not use any cleaning fluid.

- Do not use organic solvents (e.g. alcohol, ether, acetone, xylol or other dilutions) to clean lacquered parts or plastic parts!
- Disassemble optical elements (e.g. objective, eyepiece, etc) before cleaning.
- Blow away loose dust from the lens surfaces first.
- Use high quality lens cleaning towels or soft cloth and moisten it with a little bit of pure alcohol (available in a drug store). Wipe the lens surface with it.
- Pull a dust protection cover* over the microscope and store it in a closed container at a dry and mould-free place. Store the microscope and the accessories in the relevant containers when they are not used for a longer time. It is recommended to store also objectives and eyepieces in closed containers with desiccant.

(*depending on the model a dust protection cover and box can be part of the content)

NOTICE

Remember:

A well maintained microscope will keep its optical quality for years and thus maintain its value.

23 Disposal

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please make sure to dispose of your used batteries as required by law — at a local collection point or in the retail market. Disposal in domestic waste violates the Battery Directive.

Batteries that contain toxins are marked with a sign and a chemical symbol. "Cd" = cadmium, "Hg" = mercury, "Pb" = lead.

24 Technical data

	Science TFM-201	Science TFM-301	Science Infinity
Article No.	5750800	5750900	5760700
Microscope head	binocular	trinocular (incl. photo adapter)	trinocular (incl. photo adapter)
Cross table	Yes, with Nonius scale	Yes, with Nonius scale	Yes, with Nonius scale
Eyepiece(s)	1 Pair of DIN WF 10x	1 Pair of DIN WF 10x	1 Pair of DIN WF 10x
Objective(s)	4x / 10x / 40x / 100x	4x / 10x / 40x / 100x	4x / 10x / 40x / 100x
Magnification	40x / 100x / 400x / 1000x	40x / 100x / 400x / 1000x	40x / 100x / 400x / 1000x
Power supply	DC 5V / 1000mA	DC 5V / 1000mA	DC 5V / 1000mA
Illumination	LED*	LED*	LED*

*The LED illumination used in this device has a very long lifespan and do not need to be changed. LEDs can not be changed due to technical reasons!

25 Warranty

The regular guarantee period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at **www.bresser.de/warranty_terms**.

26 EC Declaration of Conformity

 Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards.

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GB IE

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail.

e-mail: service@bresseruk.com
Telephone*: +44 1342 837 098

BRESSER UK Ltd
Unit 1 starborough Farm,
Starborough Road, Nr Marsh Green,
Edenbridge, Kent TN8 5RB
Great Britain

*Number charged at local rates in the UK (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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Folux B.V.